



HEDETECH

# Odor Sampling Bag

**The Triangular Odor Bag Method<sup>1</sup>** The sensory test is conducted by at least 6 members of the panel. Each panel is given 3 bags; 1 with a sample in it and 2 without sample (odor-free air) and asked to choose the odorous bag. If the panel can tell the correct bag, the odor is then diluted and the test is continued until it becomes impossible to identify the bag with odor.



## PET odor gas sampling bag with PC stopcock valve with silicone septum for sampling the sample for the triangular odor bag method

Our odor bag come with patented valves. The materials of the valves are Polycarbonate (PC) or Polytetrafluoroethylene (PTFE) both valves have on/off function and a sample cap with septum inside for syringe sampling purpose, both valves are easy to use.

### POL Series odor bag specification

Description	POL series
Material	PET (Polyethylene terephthalate)
Appearance	Transparent
Thickness	38 $\mu$ m
Capacity	1, 3, 5, 10, 15, 20, 25, 30, 50, 80, 100 Liters
Valve type	PC Valve (Polycarbonate) -POL11 series PTFE Valve (Polytetrafluoroethylene) – POL31 series
Valve size	Inlet 6.35 mm outside diameter
Sample cap	Silicone septum for syringe sampling
Package	10 bags/box



**HEDETECH**

# Odor Sampling Bag

## Polyester odor gas sampling bag with glass tube

Concern to the triangular odor bag method. Hede Tech provide odor bag which following the method. The material made from Polyester film hot welding seal and glass tube for inlet and outlet of the sample for the panel test with 3 Liters capacity.

Description	POLOD-3L
Material	PET (Polyethylene terephthalate)
Appearance	Transparent
Thickness	38um
Capacity	3 Liters
Inlet/Outlet	10mm ID ,12mm OD Glass tube
Package	300 bags/box



### Silicone plug

- 10 mm. inside diameter
- 100 pcs/pack



### Olfactory Measurement Operator<sup>1</sup>

An olfactory measurement operator is a person in charge of management and organization of the entire series of olfactory measurement from panel selection, sampling, performance of tests and summarizing the results based on the Triangular Odor Bag Method.

### Hede Tech Co., Ltd.

6-19, 495# Changjiang Road, Xigang District,  
Dalian, Liaoning, China 116033

Tel : 86—18698631865

E-mail : sales@hedetech.com